#120

# Analysis of algae using Pyrolysis and THM

Analytical pyrolysis provides a means to analyze complex materials, even biological samples, using GCMS. Microbes, tobacco, plant fibers and foods have all been studied using this technique. Bio- fuel sources, including vegetable oils, wood, lignin and grasses have also been analyzed using Py-GC/MS, which uses thermal energy to convert large molecules like polysaccharides and proteins into smaller, volatile molecules suitable for GCMS.

Algae samples may be analyzed using a variety of thermal techniques. Figure 1 shows a sample of dried algae pyrolyzed at 600°C directly to the GC. At this temperature, the biopolymers are degraded and the pyrogram consists of characteristic frag- ments, including aromatics, nitriles and aliphatics in addition to compounds desorbed intact. Since they are a potential source of fuel, algae are fre- quently analyzed for their oil content. By pyrolysis, the fatty acids are generally fragmented, producing the normal alkanes and alkenes seen in the pyrogram.

Another approach is thermally-assisted hydrolysis/ methylation (THM) in which a reagent like tetramethyl ammonium hydroxide (TMAH) is added to the sample, which is then warmed to produce the methyl esters of the fatty acids. Figure 2 shows a sample of algae heated to 400°C in a quartz tube to which 2 µL of TMAH solution (25% in methanol) was added. The sample was allowed to stand for 5 minutes after the TMAH was added, prior to analy- sis. Under these conditions, the biopolymers are not degraded, so the resulting chromatogram is simpler, consisting almost exclusively of the fatty acid methyl esters released from the algae.

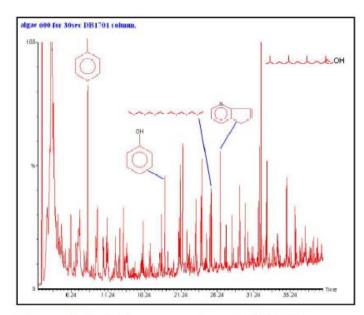


Figure 1. Dried algae pyrolyzed at 600°C.

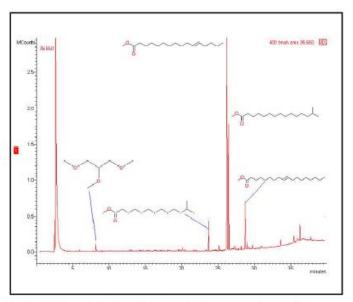


Figure 2. THM analysis of algae using TMAH.

#### **Equipment**

Samples were analyzed using a Pyroprobe 5250 Autosampler interfaced to a GC/MS. Samples were placed onto the top of a quartz spacer rod, then topped with a small plug of quartz wool to contain the TMAH solution.

## **Pyrolysis**

CDS Analytical Pyroprobe 5250 Autosampler

Pyrolysis: 600°C for 15 seconds THM: 400°C for 20 seconds

Valve oven: 300°C Transfer line: 300°C

## **GCMS**

Carrier gas: Helium Split: 50:1

Column: 25 m x 0.25 mm 5% phenyl

Oven

Initial: 40°C for 2 minutes

Ramp: 10°C/minute

Final: 300°C for 5 minutes

Range: 35 to 550

FOR MORE INFORMATION
CONCERNING THIS APPLICATION,
WE RECOMMEND THE FOLLOWING
READING:

Challinor, J. M., Review: The development and application of thermally assisted hydrolysis and methylation reactions, J. Anal. Appl. Pyrolysis, 61 (2001) 3-34.

© courtesy of CDS Analytical Inc.

#### Headquarters

JSB International Tramstraat 15 5611 CM Eindhoven T +31 (0) 40 251 47 53 F +31 (0) 40 251 47 58

Zoex Europe Tramstraat 15 5611 CM Eindhoven T +31 (0) 40 257 39 72 F +31 (0) 40 251 47 58

#### Sales and Service

Netherlands Apolloweg 2B 8239 DA Lelystad T +31 (0) 320 87 00 18 F +31 (0) 320 87 00 19

Belgium Grensstraat 7 Box 3 1831 Diegem T +32 (0) 2 721 92 11 F +32 (0) 2 720 76 22 Germany Max-Planck-Strasse 4 D-47475 Kamp-Lintfort T +49 (0) 28 42 9280 799 F +49 (0) 28 42 9732 638

UK & Ireland Cedar Court, Grove Park Business Est. White Waltham, Maidenhead Berks, SL6 3LW T +44 (0) 16 288 220 48 F +44 (0) 70 394 006 78

info@go-jsb.com www.go-jsb.com

